

QuickGuide

QuickTox Kit for QuickScan - Zearalenone

Sample Preparation



Collect and grind a representative corn sample to a 20 mesh screen size



Ground too coarse = improper extraction



Ground too fine = extract may require longer settling time

Test Procedure

(more detailed instructions in the Product Insert)




1. Add a 20 to 50 gram sub-sample to container



2. Add 2 mL/gram of 50% ethanol (example: 40 mL for 20 gram sample)



3. Shake 1 minute on mechanical shaker or by hand for 1½ - 2 minutes

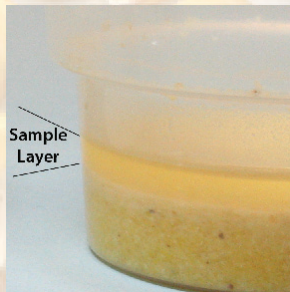

4. Allow to settle into two layers (sample is taken from top layer) and set out a vial



5. Add 100 µL Buffer to vial, then using a new pipette tip, add 100 µL sample to vial - stir well with pipette



6. Add test strip; wait 5 minutes for results

QuickScan Test

Results

(Read single strip, multiple strips, or single strip with a QuickComb—more detailed instructions in the QuickScan User Manual)

7. Remove strip from vial immediately after 5 minutes. Cut off and discard bottom pad with arrow tape. (No drying step!)

8. Place in the QuickScan carrier and slide carrier in. Click "Read Test" on Main Menu. Zearalenone results are reported between 50 ppb and 520 ppb.

9. Results Screen will appear when scanning is complete. Enter sample identification data and use buttons to save or print report.